



NeoStain ABC Kit, AP, Mouse & Rabbit

Alkaline Phosphatase labeled streptavidin-biotin detection system for broad spectrum

NB-23-00023-1

NB-23-00023-2

NB-23-00023-3

NB-23-00023-4

NB-23-00023-5

NB-23-00023-6



NeoStain ABC Kit, AP, Mouse & Rabbit

Cat# NB-23-00023-1; NB-23-00023-2; NB-23-00023-3; NB-23-00023-4; NB-23-00023-5; NB-23-00023-6

Intended Use:

Storage: 4 - 8°C

NeoStain ABC Kit, AP, Mouse & Rabbit uses biotinylated secondary antibody and Alkaline Phosphatase (AP) labeled-streptavidin to detect mouse and/or rabbit primary antibody (user-supplied) that bind to antigens in human tissue or cell preparations under light microscopy. The most commonly used specimens for this system are: frozen tissue, paraffin-embedded tissue, freshly prepared lymphocytes and fixed culture cells. Alkaline Phosphatase (AP) labeled-streptavidin and biotinylated secondary antibody amplification system has become a standard technique in immunochemical staining.

NeoStain ABC Kit, AP, Mouse & Rabbit uses human-absorbed, biotinylated, affinity-purified secondary antibody reacts with the user supplied primary antibody bound to the specific epitope of the antigen in tissue or cell. Alkaline Phosphatase (AP) labeled streptavidin then reacts with biotinylated secondary antibody to form an AP-streptavidin-biotin complex. The AP enzyme of the streptavidin complex catalyzes the substrate/chromogen such as Fast-Red, AP-Red, or BCIP/NBT to form a red (Fast-Red or AP-Red) or dark blue/purple (BCIP/NBT) color deposit at the antigen site. The antigen then can be visualized under microscope. Compared to traditional ABC methods which uses avidin, NeoStain ABC Kit, AP, Mouse & Rabbit demonstrates stronger binding strength to bind biotin and less non-specific background staining.

Higher sensitivity and lower background give this kit a higher signal-noise ratio. It also provides users cost effective method for their research. End users may choose Fast-Red, AP-Red, or BCIP/NBT chromogen depending on their preferences

Kit Components:

Reagent number:					
	1	2	3	4	
Catalog No	NeoStain ABC Kit, AP, Mouse & Rabbit	Pre-Blocking Solution	Biotinylated second antibody broad spectrum	Streptavidin-AP conjugate	Chromogen
NB-23-00023-1	no chromogen	110 ml	110 ml	110 ml	Not included
NB-23-00023-2	no chromogen	60ml	60 ml	60 ml	Not included
NB-23-00023-3	with Fast Red	8 ml	18 ml	18 ml	a. 15 Fast Redtablets b. 80ml Substrate buffer (RTU)
NB-23-00023-4	with Fast Red	6 ml	6 ml	6 ml	a. 6 Fast Red tablets b. 35ml Substrate buffer (RTU)
NB-23-00023-5	with Permanent Red	18 ml	18 ml	18 ml	a. 1.5ml AP-Red+ Enhancer (40x) b. 1.5ml AP-Red+ Solution (40x) c. 6ml AP-Red+ Substrate (20x)
NB-23-00023-6	with Permanent Red	6 ml	6 ml	6 ml	a. 1ml AP-Red+ Enhancer (40x) b. 1ml AP-Red+ Solution (40x) c. 4ml AP-Red+ Substrate (20x)

Recommended Protocol:

1. Fixation: To ensure the quality of the staining and obtain reproducible performance, user needs to supply appropriately fixed tissue and well prepared slides.
2. Tissue need to be adhered to the slide tightly to avoid tissue falling off.
3. Paraffin embedded section must be deparaffinized with xylene and rehydrated with a graded series of ethanol before staining.
4. Cell smear samples should be made as much monolayer as possible to obtain satisfactory results.
5. Three control slides will aid the interpretation of the result: positive tissue control, reagent control (slide treated with Isotype control reagent), and negative control.
6. Start staining procedures: DO NOT let specimen or tissue dry from this point on.

Reagent	Staining Procedures	Incubation Time (Min.)
1. HIER Pretreatment: refer to antibody spec. sheet	a. Heat Induced Epitope Retrieval (HIER) may be required for primary antibody suggested by vendor. b. Wash with PBS 3 times for 2 minutes each time.	
2. Reagent 1: Pre-blocking Solution	a. Add 2 drops or enough of volume Pre-blocking Solution to completely cover the tissue section and Incubate for 10 min. b. Blot off solution. DO NOT RINSE.	10 min.
3. Primary antibody: Supplied by user. Investigator needs to optimize dilution and incubation time	a. Apply 2 drops or enough volume of Primary antibody to cover the tissue section completely. Incubate in moist chamber for 30-60 min. b. Rinse with PBS 3 times for 2 minutes each time.	30-60 min.
4. Reagent 2: Ready-to-use Broad Secondary antibody	a. Apply 2 drops or enough volume of secondary antibody to cover the tissue section completely and incubate for 10 min. b. Rinse with PBS 3 times for 2 minutes each time	10 min.
5. Reagent 3: Ready-to-use AP-Streptavidin	a. Apply 2 drops or enough volume of AP-Streptavidin to cover the tissue section completely and incubate for 10 min. b. Rinse with PBS 3 times for 2 minutes each time. c. Rinse with tap water.	10 min.
6. Reagent 4: Chromogen: Fast-Red, or AP-red, or BCIP.NBT	Refer to manufacture data sheet if chromogen is supplied by user. Recommended protocol for chromogen using our kit: 1. Fast Red : a. Dissolve one Fast Red tablet into one 5ml substrate buffer. Vortex until tablet is dissolved. It usually takes 20 minutes to dissolve completely. b. Chromogen must be used within 1 hour. c. Apply 100ul or more Fast-Red solution to completely cover the tissue section and incubate 10 minutes at room temperature. d. After proper color development, wash with distill water for 2 minutes, 3 times e. DO NOT Dehydrate tissue after staining. Fast-Red is alcohol soluble. 2. AP-Red+ (40x):	10 min.

	<p>a. Add 1 drop (50ul) of AP-Red+ Enhancer and 1 drop (50ul) of AP-Red+ Solution to a test tube. Mix well and set at room temperature for about 5 minutes.</p> <p>b. Add 2ml of distilled water to the mixture. Mix well.</p> <p>c. Add 4 drops (200ul) of AP-Red+ Substrate to the mixture and mix well.</p> <p>d. Completely cover the tissue section with the mixture and incubate for 5-15 minutes.</p> <p>e. After proper color development, wash with distill water 2 minutes, 3 times. APRed+ is soluble in organic solvent. Do not dehydrate.</p> <p>3. BCIP/NBT : order separately, Cat. No. NB-23-00144-1 or NB-23-00144-2</p> <p>a. Add two drops (about 100ul) of Ready-to-use BCIP/NBT to cover the tissue section for 5-10 minutes. Monitor the color development under a microscope. b. Rinse with distill water for 2 minutes, 3 times</p>	
<p>7. Hematoxylin: Supplied by user</p>	<p>a. Counterstain with 2 drops or enough volume to cover tissue completely and wait about 10-20 seconds.</p> <p>b. Rinse thoroughly under tap water for 1-2 min.</p> <p>c. Put slides in PBS until show blue color (about 30-60 seconds) d. Rinse well in distilled water</p>	
<p>8. Mounting media: Supplied by user</p>	<p>Follow the manufacturer's data sheet procedure for mounting. Recommended product:</p> <p>1. NeoMount AQ : Cat. No. NB-23-00155-3 (18ml) for Aqueous solution, use with AEC and AP-Red</p> <p>2. NeoMount Perm: Cat. No. NB-23-00156 (18ml), for Permanent mount for DAB,BCIP/NBT</p> <p>3. NeoMount Universal: Cat.No. NB-23-00157-2 (18ml), or NB-23-00157-1 (100ml), for Water Based, universal</p>	

Protocol Notes:

1. The fixation, tissue slide thickness, antigen retrieval and primary antibody dilution and incubation time effect results significantly. Investigator needs to consider all factors and determine optimal conditions when interpret the result.
2. Tissue staining is dependent upon the proper handling and processing of tissues prior to staining. Improper tissue preparation may lead to false negative results or inconsistent results.
3. Do not mix reagents from different lot.
4. Do not allow the slides to dry at any time during staining

Related Products:

Product	Catalog No	Size	Product	Catalog No	Size
Mouse , no chromogen	NB-23-00024-1	110 ml	Fast Red Kit	NB-23-00142	12 Tab + 60ml
Mouse , with Fast Red	NB-23-00024-2 NB-23-00024-3	18ml / 6ml	AP-Red concentrated (40x) Kit for 2000 slides	NB-23-00143	8ml
Mouse , with Permanent Red	NB-23-00024-4 NB-23-00024-5	18ml / 6ml	BCIP/NBT (RTU)	NB-23-00144	100ml / 18ml
Rabbit , no chromogen	NB-23-00025-1	110ml	NeoMount AQ (Aqueous solution, use with AEC and AP-Red)	NB-23-00155-3	18ml
Rabbit , with Fast Red	NB-23-00025-2 NB-23-00025-3	18ml / 6ml	NeoMount Perm (Permanent mount for DAB,BCIP/NBT)	NB-23-00156	18ml
Rabbit , with Permanent Red	NB-23-00025-4 NB-23-00025-5	18ml / 6ml	NeoMount Universal (Water Based, universal) Kit	NB-23-00157-	100ml / 18ml
Streptavidin-AP (RTU)	NB-23-00027-1 NB-23-00027-2	100ml / 18ml			

Precautions:

Handle all specimens as potential infectious materials, wear gloves and protection cloth.

Remarks:

For research use only