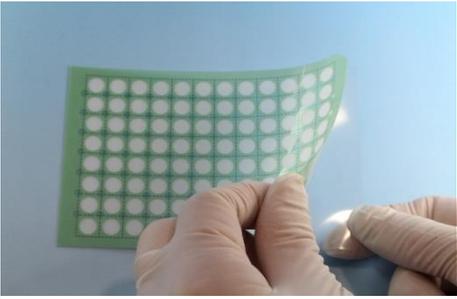
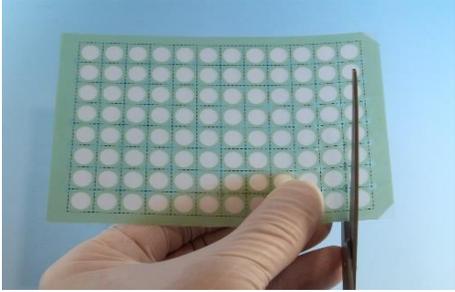
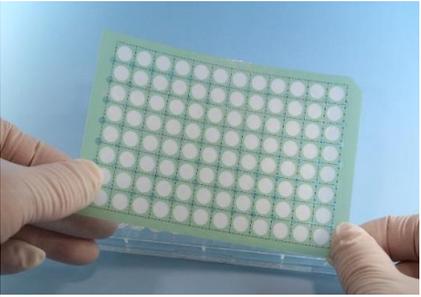
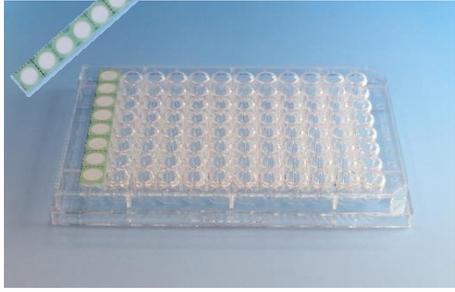
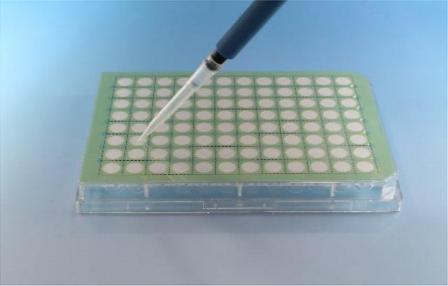


Directions of Microplate Mesh Seal for 96-well plate

	96 well use		8 well use
①	 <p>Peel off the liner on the back side.</p>	①	 <p>Tear off along the perforation by scissors. Possible to tear off just 1 well.</p>
②	 <p>Place the mesh seal on the microplate along the well position. Possible to replace many times.</p>	②	 <p>Place the torn off one on the microplate after peeling off the liner.</p>
③	 <p>Put drops of specimen by pipette from an angle.</p>	③	 <p>Put drops of specimen by pipette from an angle.</p>