

Foxx Life Sciences' Heavy Duty (HD) Solvent Waste System is designed to be an affordable solution for liquid waste disposal. Each system is manufactured to safely collect waste from one or more HPLC systems while the carbon exhaust filter improves air quality by capturing potentially harmful vapors emitted from solvents. Each system includes a 2.5L, 4L, 5L, 10L, or 20L HDPE container, two 83B **VersaCap®** (one open top for waste collection and one closed for safe transport), various ports for waste tubing, 1 port for a carbon exhaust filter, 1 carbon exhaust filter, and plugs for unused ports.



The HD Solvent Waste System manifold attaches to the **EZgrip™** Carboy allowing for safe waste collection.

- 1** Carbon exhaust filter* traps solvent vapors
- 2** Tube fittings secure waste tubes to carboy providing a closed system that prevents vapors from escaping
- 3** 90° adapter included for maximum storage flexibility
- 4** **VersaCap®** technology eliminates the need to remove tubing from open cap to install closed cap
- 5** Plugs included for unused ports
- 6** Replacement sulfur-cured EPDM gasket for high chemical compatibility against aggressive organic solvents
- 7** Closed 83B **VersaCap®** included for safe transport
- 8** High Density Polyethylene **EZGrip™** carboy features advanced ergonomic design and superior chemical compatibility

*Depending on application, carbon exhaust filter life cycle may vary



Distributed by:

CliniSciences Group ISO 13485:2016




	1/8" OD PORTS	1/4" OD PORTS	1/4" or 3/8" HB PORT	PART NUMBER
2.5L	6	-	-	332-1412-OEM
	6	-	1	332-1418-OEM
	4	3	-	332-1413-OEM
	4	4	-	332-1421-OEM
	4	3	1	332-1419-OEM
4L	6	-	-	332-9412-OEM
	6	-	1	332-9418-OEM
	4	3	-	332-9413-OEM
	4	4	-	332-9421-OEM
	4	3	1	332-9419-OEM
5L	6	-	-	332-2412-OEM
	6	-	1	332-2418-OEM
	4	3	-	332-2413-OEM
	4	4	-	332-2421-OEM
	4	3	1	332-2419-OEM
10L	6	-	-	332-3412-OEM
	6	-	1	332-3418-OEM
	4	3	-	332-3413-OEM
	4	4	-	332-3421-OEM
	4	3	1	332-3419-OEM
20L	6	-	-	332-4412-OEM
	6	-	1	332-4418-OEM
	4	3	-	332-4413-OEM
	4	4	-	332-4421-OEM
	4	3	1	332-4419-OEM
20L ROUND	6	-	-	332-D412-OEM
	6	-	1	332-D418-OEM
	4	3	-	332-D413-OEM
	4	4	-	332-D421-OEM
	4	3	1	332-D419-OEM

* 1/16" OD Adapters are also available for the HD line as cap only assemblies.


Solvent Waste System Sizing Guide

Due to all the possibilities with configuring your EZwaste Solvent Waste System, we have provided the following images to help identify the fitting required for your laboratory setup. All measurements are approximate.

① For measuring OD, place the tube next to each of the images to determine the correct size of the tubing.

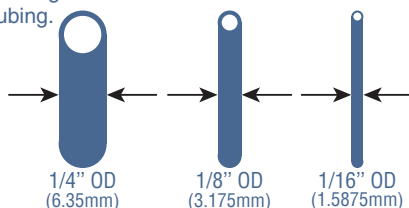

② For measuring ID, place end of tube flat on the piece of paper and determine if the dark area is contained by the tubing.

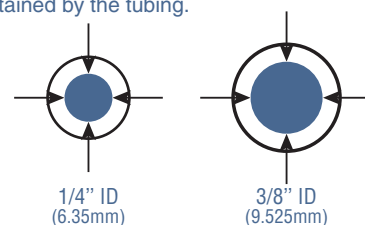

CliniSciences Group

Austria



Company: CliniSciences GmbH
Address: Sternwartestrasse 76, A-1180
Wien - Austria
Telephone: +43 720 115 580
Fax: +43 720 115 577
Email: oesterreich@clinisciences.com
Web: <https://www.clinisciences.com>

Belgium



Company: CliniSciences S.R.L
Address: Avenue Stalingrad 52, 1000
Brussels - Belgium
Telephone: +32 2 31 50 800
Fax: +32 2 31 50 801
Email: belgium@clinisciences.com
Web: <https://www.clinisciences.com>

Denmark



Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: denmark@clinisciences.com
Web: <https://www.clinisciences.com>

Finland



Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: suomi@clinisciences.com
Web: <https://www.clinisciences.com>

France



Company: CliniSciences S.A.S
Address: 74 Rue des Suisses, 92000
Nanterre- France
Telephone: +33 9 77 40 09 09
Fax: +33 9 77 40 10 11
Email: info@clinisciences.com
Web: <https://www.clinisciences.com>

Germany



Company: Biotrend Chemikalien GmbH
Address: Wilhelm-Mauser-Str. 41-43,
50827 Köln - Germany
Telephone: +49 221 9498 320
Fax: +49 221 9498 325
Email: info@biotrend.com
Web: <https://www.biotrend.com>

Iceland



Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: island@clinisciences.com
Web: <https://www.clinisciences.com>

Ireland



Company: CliniSciences Limited
Address: Ground Floor, 71 lower Baggot street
Dublin D02 P593 - Ireland
Telephone: +353 1 6971 146
Fax: +353 1 6971 147
Email: ireland@clinisciences.com
Web: <https://www.clinisciences.com>

Italy



Company: CliniSciences S.r.l
Address: Via Maremmana inferiore 378
Roma 00012 Guidonia Montecelio - Italy
Telephone: +39 06 94 80 56 71
Fax: +39 06 94 80 00 21
Email: italia@clinisciences.com
Web: <https://www.clinisciences.com>

Netherlands



Company: CliniSciences B.V.
Address: Kraijenhoffstraat 137A,
1018RG Amsterdam, - Netherlands
Telephone: +31 85 2082 351
Fax: +31 85 2082 353
Email: nederland@clinisciences.com
Web: <https://www.clinisciences.com>

Norway



Company: CliniSciences AS
Address: c/o MerVerdi Munkerudtunet 10
1164 Oslo - Norway
Telephone: +47 21 988 882
Email: norge@clinisciences.com
Web: <https://www.clinisciences.com>

Poland



Company: CliniSciences sp.Z.o.o.
Address: ul. Rotmistrza Witolda Pileckiego 67
lok. 200 - 02-781 Warszawa -Poland
Telephone: +48 22 307 0535
Fax: +48 22 307 0532
Email: polska@clinisciences.com
Web: <https://www.clinisciences.com>

Portugal



Company: Quimigen Unipessoal LDA
Address: Rua Almada Negreiros, Lote 5, Loja 14,
2615-275 Alverca Do Ribatejo - Portugal
Telephone: +351 30 8808 050
Fax: +351 30 8808 052
Email: info@quimigen.com
Web: <https://www.quimigen.pt>

Spain



Company: CliniSciences Lab Solutions
Address: C/ Hermanos del Moral 13
(Bajo E), 28019, Madrid - Spain
Telephone: +34 916 750 700
Fax: +34 91 269 40 74
Email: espana@clinisciences.com
Web: <https://www.clinisciences.com>

Sweden



Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: sverige@clinisciences.com
Web: <https://www.clinisciences.com>

Switzerland



Company: CliniSciences AG
Address: Fracht Ost Flughafen Kloten
CH-8058 Zürich - Switzerland
Telephone: +41 (044) 805 76 81
Fax: +41 (044) 805 76 75
Email: switzerland@clinisciences.com
Web: <https://www.clinisciences.com>

UK



Company: CliniSciences Limited
Address: 11 Progress Business center, Whittle
Parkway, SL1 6DQ Slough- United Kingdom
Telephone: +44 (0)1753 866 511
or +44 (0) 330 684 0982
Fax: +44 (0)1753 208 899
Email: uk@clinisciences.com
Web: <https://www.clinisciences.com>

USA



Company: Biotrend Chemicals LLC
Address: c/o Carr Riggs Ingram,
500 Grand Boulevard, Suite 210 Miramar
Beach, FL 32550- USA
Telephone: +1 850 650 7790
Fax: +1 850 650 4383
Email: info@biotrend-usa.com
Web: <https://www.biotrend-usa.com>